

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922) 49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Ижевск (3412)26-03-58
Иваново (4932)77-34-06
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Киргизия (996)312-96-26-47

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Россия (495)268-04-70


Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сыктывкар (8212)25-95-17
Сургут (3462)77-98-35
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Казахстан (772)734-952-31

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Улан-Удэ (3012)59-97-51
Ульяновск (8422)24-23-59
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

<https://integra.nt-rt.ru> || iag@nt-rt.ru

Наконечники автоматизированные для пипеток GRIPTIPS

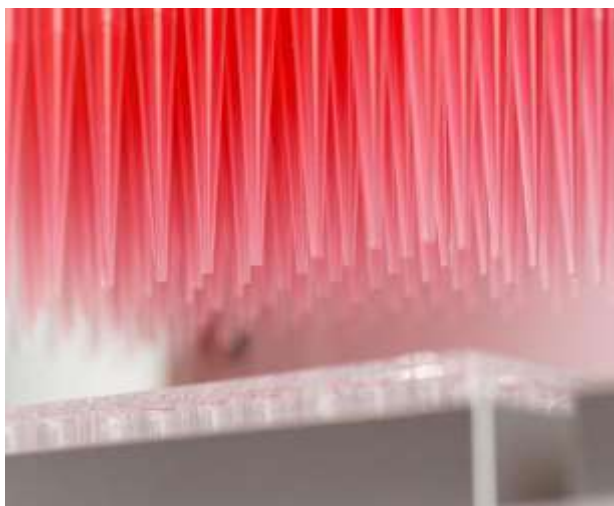

Perfecting the connection between pipetting head and tips

Do your pipette tips ever loosen, leak or fall off? This is a common issue in laboratories, caused by the use of universal pipette tips. Losing a tip during assay preparation can ruin the entire assay; a reliable seal is essential for accurate pipetting.

INTEGRA has redesigned the connection between the pipette tip and the tip fitting to resolve this issue. GRIPTIP pipette tips perfectly match the multi-lobe tip fitting, snapping firmly on during loading. This guarantees a perfect seal on every tip, preventing them from loosening, leaking or completely falling off. All tips are precisely aligned horizontally, enabling accurate touch-offs, even when pipetting with 384 tips.

GRIPTIPS suitable for our benchtop pipetting platforms undergo additional quality control procedures – such as a secondary straightness test – to guarantee perfect targeting of wells for automatic pipetting or simultaneous pipetting with up to 384

tips. The tip racks are also reinforced to withstand the forces required to load 384 tips.

How it works

Precise and consistent tip seal ensures that all tips on VIAFLO 96, VIAFLO 384, and ASSIST PLUS are precisely at an identical height for reliable touch offs and consistent pipetting results. An additional tip straightness inspection process ensures all pipette tips access wells even when dispensing to a 384 well plate.

• **GRIPTIP Pipette tips which never leak or fall off**



Losing a tip during an assay preparation can ruin the entire assay and a reliable sealing is of key importance for accurate pipetting.

INTEGRA has redesigned the connection between pipette tip and tip fitting to resolve this issue. The unique GRIPTIP pipette tips are comfortably loaded and ejected without any physical effort due to a power assisted mechanism activated by a push button. The loading procedure for all benchtop pipetting platforms also guarantees that GRIPTIPS are not touched by the user and remain sterile.

-
- 1) LOBES
 - 2) O-RING SEAL
 - 3) RIM SNAPS OVER THE LOBES
 - 4) SHOULDER
-

Extended length design



INTEGRA's GRIPTIPS LONG feature an extended length design enabling easy access into labware types such as tubes and deep well blocks.

The extended length design also reduces PCR cross contamination as the elongated tip prevents the pipette shaft from contacting the wall of the PCR tubes.

Low retention GRIPTIP pipette tips



Viscous samples, detergents and other low surface tension liquids tend to interact with standard pipette tips forming a thin liquid film

on the pipette tip inner wall. This effect leads to pipetting inaccuracy and inconsistency as well as loss of precious reagents. INTEGRA uses a unique polypropylene blend to mold Low Retention GRIPTIPS with heightened hydrophobic properties. Using Low Retention GRIPTIPS prevents low surface tension samples from spreading out and “wetting” the inner wall of the tips, allowing them to bead-up for a maximum liquid recovery.

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922) 49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Ижевск (3412)26-03-58
Иваново (4932)77-34-06
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Киргизия (996)312-96-26-47

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Россия (495)268-04-70

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сыктывкар (8212)25-95-17
Сургут (3462)77-98-35
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Казахстан (772)734-952-31

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Улан-Удэ (3012)59-97-51
Ульяновск (8422)24-23-59
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

<https://integra.nt-rt.ru> || iag@nt-rt.ru